Dino-Lite

Digital Microscope

The Industry Standard

Dino-Lite

User Manual

EN	Manual	PT	Manual do Utilizador
FR	Manuel d'utilisation	RU	Руководство по ЭКСПЛУАТАЦИИ
DE	Bedienungsanleitung	SE	Användarhandbok
ES	Manual de usuario	PL	Instrukcja obsługi
IT	Manuale utente	TR	Kullanım Kılavuzu
NL	Handleiding		

www.dino-lite.eu

Redefine the microscope

Dino-Lite hardware and software user manual

ENGLISH	3
FRANCAIS	15
DEUTSCH	27
ESPAÑOL	39
ITALIANO	51
NEDERLANDS	63
PORTUGUÊS	75
РУССКИЙ	88
SVENSKA	102
TÜRKÇE	114
POLSKI	126

English

Thank you for purchasing a Dino-Lite or Dino-Eye digital microscope. The DinoCapture software is designed to give you the best possible digital microscopy experience by the inventors of the handheld digital microscope. The Dino-Capture software runs on computers with a Windows XP, Windows Vista or Windows 7 /8 operating system. For Apple Macintosh OS, please refer to the DinoXcope software that can be downloaded from www.dino-lite.eu.

⚠ Important safety information

- Avoid touching the lens with finger to protect the product from electrostatic damage.
- Do not drop
- Keep dry

Index

Installation	4
Software interface (DinoCapture / Windows)	6
Software interface (DinoXcope / Mac)	11
Cleaning & Maintenance, Warranty, Support	13

Software installation

Important notice: DO NOT connect the USB cable of the Dino-Lite or Dino-Eye to the PC before installing the software

1. Use the CD delivered with your Dino-Lite to install the DinoCapture software and drivers. Alternatively, download and run the latest version of the software from the support section of the website: www.dino-lite.eu.
2. Click 'Next' and the Installshield wizard will start. (An 'Open File-security warning' may appear on some systems. Select 'Run' or 'YES'). Choose the language you want for the DinoCapture 2.0 interface.
3. Read the User License Agreement. If you agree, press 'Yes' to continue or press 'no' to stop the installation.
4. Select a destination folder for the DinoCapture software. When done, press 'Next'. Click 'Install' to start installing the software. If the Windows security warning message appears, click 'Install this driver software anyway'.
5. When the installation is complete, selecting 'Finish' completes the software installation.
6. The DinoCapture software has an auto-update feature that will check for software updates when you start DinoCapture.
7. A full manual can be found in the help function of DinoCapture, as a pdf on the CD or on the website.

Hardware installation

1. After full installation of the DinoCapture software and driver package, connect the Dino-Lite to one of the USB ports of your computer.
2. Please use a USB 2.0 port that is fully powered. Some USB ports on portable computers do not supply sufficient power.
3. The driver will be automatically installed. Please WAIT until the notification appears: 'Device driver software installed successfully'.
4. Now start DinoCapture 2.0 by double clicking on the desktop icon.

5. The LED lights should go on and an image should appear in DinoCapture. If this is not the case, refer to the frequently-asked-questions (FAQ) on www.dino-lite.eu.

Hardware features

- a. At the center of the device, the adjustable dial is used to set the focus. The focus of the image depends on the distance to the object. Once you have focused on the object, you can read the magnification rate achieved from the number next to the Δ symbol.
- b. The models in the AM/AD 411X series feature a magnification lock. The magnification lock is particularly useful for repetitive inspection at a given level of magnification.
- c. Dino-Lite models with the letter T in their product name have a MicroTouch function at the cable end of the device. Touching this sensor will capture the current image or start/stop video recording (for usb-devices). On the Dino-Lite high speed real time models the Microtouch sensor can be used to switch the LEDs on/off (all models) or freeze the image (AM5116 models only).
- d. Models with the letter Z in their product name have a polarization function, that can be controlled by turning the adjustable cap.
- e. Models with the letter L operate at a longer working distance. These models only achieve focus at some distance from the object.
- f. The AD models have exchangeable front covers/caps. Align the red lines on cap and frontend of the Dino-Lite to remove or place the cap, then turn the cap 180 degrees. Some caps are designed to be clicked onto the microscope. In this case, there is no red line on the cap.
- g. The DinoEye models are designed as a replacement of the existing ocular of a traditional microscope. The U model is intended to be placed over an existing ocular and the C model is intended to be connected to a C-mount adapter on a suitable microscope or optical device.

Software interface (Windows)

There are five key sections in the DinoCapture software:

1. Menu bar
2. Image tool list
3. Preview window management bar
4. Tool bar
5. Preview window tools

Menu bar

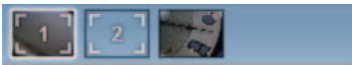
	Folder – create folders or access existing folders.
	Files – manage files, such as open, copy, save, print or delete as well as send as e-mail or create slide show.
	Settings – customize the way the software operates such as showing or hiding elements on the screen, the properties for measurements, the way the microtouch or foot pedal (if attached) work, connecting over IP and the auto update function.
	Help – access the full user guide on screen, or read DinoCapture 2.0 information or the license agreement.

Image list tools

	Open – open a picture selected from the image thumbnail gallery below the image list tool bar.
	Copy – copy a selected picture to the windows clipboard.
	Save as – save selected picture(s) to the desired file format or folder.
	E-mail – attach the selected picture(s) to a new message opened in your default e-mail client.
	Print – print the selected picture.
	Slideshow – run a slideshow of the selected pictures.
	Delete – clean up the image thumbnail gallery by deleting the selected picture(s).

Preview window management bar

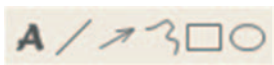
This window enables us to switch between windows if two or more Dino-Lite's are connected at the same time or multiple pictures are opened.


Tool bar

The tool bar is divided into three parts:

Draw tools - The draw tool set allows you to write text and draw on the images. The bar is located below the preview window management bar.

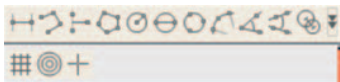Measurement tools – Many different measurement types can be done with the measurement tools. The double flash icon gives you access to the grid and ruler settings.

Notice: The measurement and calibration feature is only available for certain models. For the DinoEye models, it is recommended to make calibrated measurements only. Please contact a local sales representative for further information.

Line	Left click and drag to the desired length, and click again to finish.
Continuous line	Click and drag to form one section of distance, click again to start another section. Continue until the total desired distance is measured. Double click to finish.
Point to line	Create a line to represent the base by clicking once to start, drag, and then click again to set the endpoint. Branch off from the base line to start measuring the line that is 90 degrees from the base line and a final click to finish the measurement.

Polygon	Click and drag to form the desired length, and click again to start the next section. When finished, double click to finish the polygon measurement.
Radius	Click and extend out to the desired radius.
Diameter	Click and extend out to the desired diameter.
Three point circle	Click on any three points on the circle you wish to measure.
Three point arc	Click on three consecutive points on an arc to measure.
Three point angle	Start at the pivot point and extend out to start measuring an angle.
Four points angle	Select two points from one line segment and another two points from another line segment to complete the angle measurement.
Center distance	After drawing at least two circles, select the icon and the mouse pointer will change to a pointing finger for selectable circles. Click on each circle and the software will measure the distance.
Gridlines	The pitch will match the magnification inputted.
Circle grid	The pitch will match the magnification inputted.
Crosshair on camera	The crosshair XY position can be moved when you select it. The cursor position is the location of the mouse pointer and location 0,0 is the centre coordinates of the crosshair.
Scale crosshair	The scale crosshair increments can be compared to the scale on the bottom left corner. The centre of the crosshair can be moved. (Magnification value needs to be entered for this feature to work).

Text and line tools

Line format	Select line color, style and width
Font	Select font, font style, size and color

Measurement options

Measurement properties	Organize and show the status of all the measurements as well as the ability to adjust how the results are shown on the image.
Magnifier	Digitally enlarge the area around your mouse pointer for more details and accurate measurements.
Magnification	Input the magnification in the blue box. The magnification can be read from the dial on the microscope.
Units	Select the unit of measurement, inch, mil, mm or um.
Calibration menu	Select calibration, create new calibration profile or open the calibration folder.
Barcode Reader	Click on the icon to enable and disable barcode detection

Preview window tools

	Microtouch - this is activated when you touch the Microtouch sensor. Activate or deactivate the Microtouch. Some Dino-Lite models do not have this feature.
	Auto Exposure - This allows you to change the exposure or turn off auto exposure. Sliding the bar to the right increases the exposure time and vice versa.
	LED control - lets you turn ON/OFF or switch LED's on the Dino-Lite (most models).

	Setting - The setting window allows you to control the camera settings.
	Full screen - Lets you see the live video or picture in full screen. To exit full screen mode, click anywhere on the screen or press "ESC".
	Close window - Allows you to close the current window
Snapshot	Take a picture.
Video recorder	Start/stop recording of a video
Time-Lapsed Video	Create a series of images or videos taken at a regular time interval.
	Choose the resolution for the image.
	Click on the gray bar on the bottom of the preview window and the annotation box will expand. Write text within this box and press ENTER to start a new row. When done, left click anywhere in the preview window to save the annotation and exit the annotation box.

Software interface (DinoXcope for Mac)

There are three main sections in the DinoXcope program window:

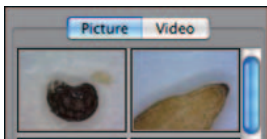
1. Action bar

- a. Make picture
 - If you want to make measurements on the picture later, open the picture by double clicking and add the input magnification that you read from the focus dial.


- b. Make video
 - Choose frame rate, recording time, quality and compression.
 - To stop recording manually, use the ESC key.
- c. Make time lapse video
 - Choose capture interval, number of frames and playback frame rate, as well as quality and compression.
- d. Switch LED's on/off
- e. Go to full screen mode, to end full screen mode use the ESC key

2. Image list

- a. Choose between pictures or video's
 - b. Look at an image with mouse-over, or double-click to open in new window
- ## 3. Live image
- a. Use the action bar for the action that you want to perform.
 - b. Use the Controls menu (the menu bar is on top of the screen) to change settings, use (digital) zoom, freeze the image, change the resolution or change the functioning of the Microtouch button.
 - c. Use the Live measurement function (in the Tools menu) to open a special selection window for drawing and measurement tools. After selection, click okay and use the tool on the live image.

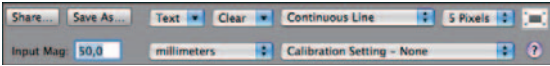4. Open image

- a. If you open a stored image, you will see a menu bar above it, that allows you to choose saving, drawing and measurement tools.

Cleaning and maintenance

Avoid touching the lens with any substance. Clean the lens periodically with compressed air. Clean the body of the microscope regularly with a soft cloth dampened with disinfectant. Clean the detachable caps with a 50%-70% alcohol solution, do not use diethyl ether, 100% alcohol or alcohol gel.

Warranty

The DinoCapture software is provided for use with a Dino-Lite or DinoEye digital microscope supplied by AnMo Electronics or one of its distributors or resellers. The use of the DinoCapture software is guided by the End User License Agreement.

The Dino-Lite product is delivered with a two year warranty from the date of purchase by the end customer. Please note that accessories (i.e. detachable caps/extensions) are not covered under warranty. For warranty issues please contact the reseller or store where you bought the product.

Support

If you have any problem or issue with your Dino-Lite or the DinoCapture software, please contact your reseller or check the Dino-Lite Europe support pages at www.dino-lite.eu. An extensive user manual for DinoCapture can be found on the website.

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e-mail: info@dino-lite.eu

web: www.dino-lite.eu

Version 2013Q3 © AnMo/IDCP/Dino-Lite Europe.

No part of this publication may be reproduced or distributed unless for personal use in conjunction with usage of Dino-Lite digital microscopes.

Patent info:

US Pat. No. 7.102.817.B1

Germany Pat. No. 20 2006 001 409.8

Japan Pat. No. 3123176

Taiwan Pat. No. M302031

AnMo Electronics Corporation and its authorized representative in the EU, IDCP BV, declare that the apparatus Dino-Lite and Dinoeye comply with the essential requirements and other relevant provisions of Directive 1999/5/EC. The original Declaration of Conformity can be found at www.dino-lite.eu/doc/dino-lite.pdf

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Version 2013Q3 Copyright © AnMo/IDCP/Dino-Lite Europe
Version 2013Q3 Copyright © 2013 IDCP BV
Dino-Lite is © of AnMo Electronics Corporation